# **ExAO-NG**

Capteur

# **ExAO-NG**

*Sensor*

Réf :  
**482 044**

Français – p 1

English – p 4

Version : 3106

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**0 825 563 563 \***

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## 1 General

This sensor, connected to CO<sub>2</sub> probe ref. 453 076, measures CO<sub>2</sub> in air or water.

## 2 Use

This sensor can be used with any console in the Foxy range.

### 2.1 Connection to the Foxy console

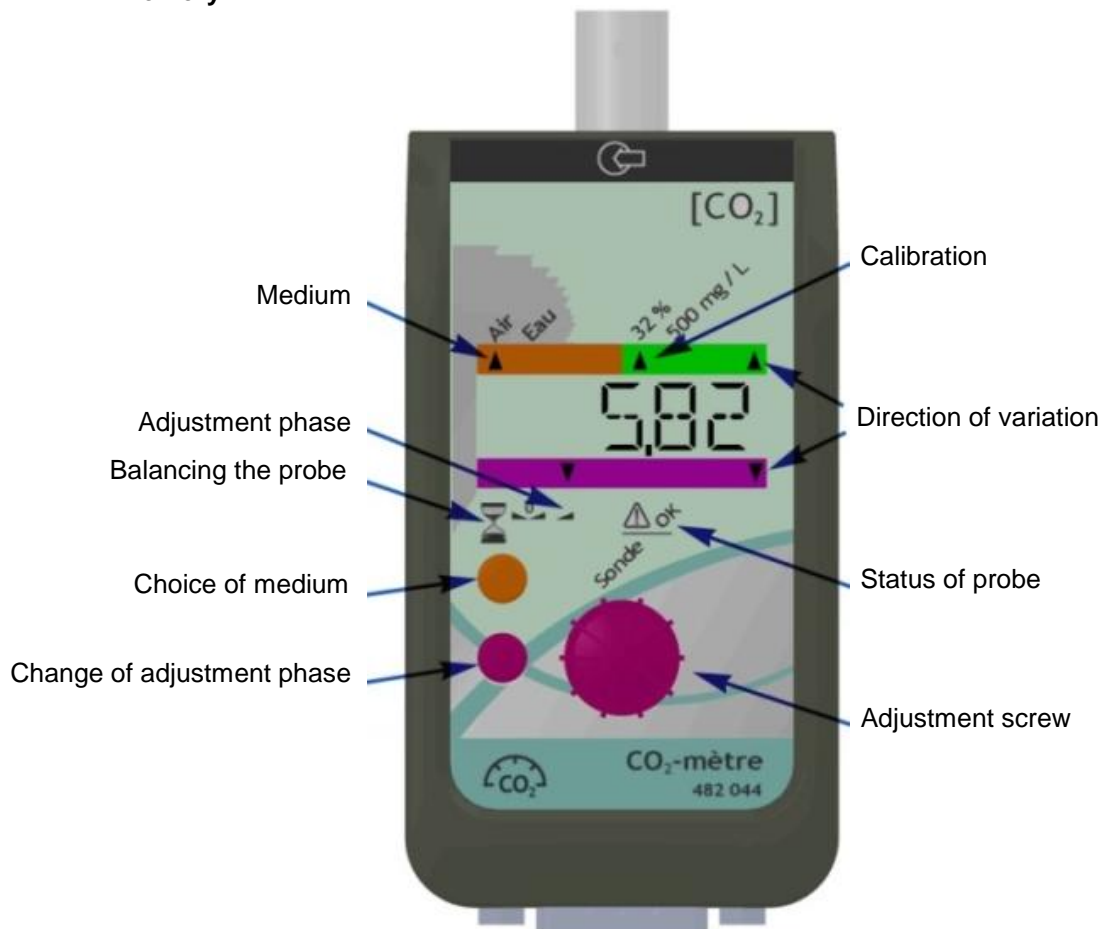
- Place the sensor opposite the console connector
- Slide the connector following the console guide
- Push home until it clips in.

### 2.2 Use in stand-alone or connected mode

#### Calibration:

#### Important remarks:

For accurate quantitative measurements, the probe must be calibrated but this is not essential every time it is used provided the probe/sensor combination remains unchanged: the sensor stores the settings in its memory.

## 2.2.1 In stand-alone mode

All the functions of a multi-meter are available. The instantaneous value of the CO<sub>2</sub> content is visible on the display.

A triangle on the display indicates the calibration and medium selected.

Adjustment:

- Wait a few minutes until the electronics are in balance: the triangle above the egg timer flashes during this phase, then stops.
- Select the medium "air" by pressing the medium selection button.
- Hold down the violet button. The triangle flashes opposite the pictogram .
- Set zero using the adjustment button, with the probe in free air.
- Briefly press the violet button. The triangle flashes opposite the "slope" pictogram .
- Place the probe in the calibration solution (see instructions for CO<sub>2</sub> probe ref. 453 076); the arrow of direction of variation at the top flashes, indicating an increase in the CO<sub>2</sub> content (N.B. the displayed value does not change).
- When the direction of variation arrows flash alternately, the probe is in balance; simply turn the adjustment button to display 5.82 (%) then press the adjustment button a final time.
- If adjustment is satisfactory, the triangle lights up opposite the "ok" pictogram.
- If the probe malfunctions, the triangle is located opposite the pictogram "HS" (malfunction).

*Remark: The fact the "malfunction" message flashes does not necessarily signify that the probe is not working. This message frequently indicates a malfunction which must be diagnosed before starting acquisition of the reading. The malfunction may result from multiple causes: pierced membrane, inadequate quantity of electrolyte, contaminated electrode, wrong electrolyte (mix up of electrolyte for the O<sub>2</sub> probe and the CO<sub>2</sub> probe, mix up between electrolyte for the CO<sub>2</sub> probe and the acid calibration solution) etc.*

## 2.2.2 In connected mode

Once the console is connected to the computer, the sensor display goes out. The values measured are visible only on the computer and the buttons become inactive. The adjustment phase is identical as in the method described above but the instructions are displayed on the computer screen.

## 3 Additional equipment

The sensor used is the air/water CO<sub>2</sub> probe ref. 453 076.

## 4 Technical characteristics

- Calibration: air 0-32%, water 0-500 mg/L
- Accuracy: 1%
- Response time at 90% : ≤ 2 min

## 5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
468, Rue Jacques Monod  
CS 21900  
27019 EVREUX CEDEX France

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

### Free service \*

**+ 33 (0)2 32 29 40 50\*\***

\* Call cost not included

\*\* Only for call from foreign countries