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**Tooxy**  
*Sensor*

Ref :  
482 204

Français – p 1

English – p 4

Version : 0112

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***Fast thermometer sensor***

## 1 Généralités

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## 1 General

This sensor is used to make temperature measurements within the  $-50^{\circ}\text{C}$  to  $+150^{\circ}\text{C}$  range.

Measurements are made using a low inertia Pt 1000 probe.

## 2 Use

This sensor operates with a console in standalone mode (without computer), or in connected mode (with computer) after being connected to the console.

### 2.1 Connection to the console

The sensor is connected using the connector on the back of the housing:

- Bring the sensor in line with the selected measurement channel
- Insert the sensor into the console guide system
- Follow the console guide system.
- Push all the way in

## 2.2 Type of measurement

The sensor measures the temperature on a scale from -50 to +150°C.

## 2.3 Operation in standalone mode

### c. Start up

- Connect the sensor
- Switch the console on

### d. Display

An instantaneous temperature value is displayed directly on the display screen present on the housing.

## 2.4 In connected mode

### c) Start up

- Connect the sensor to the console
- Switch the console on
- Connect the console to the computer

### d) Display

Measurements are displayed directly on the computer screen, either during an acquisition or using the “display” tab

## 3 Technical characteristics

Measurement range	-50/+150°C
Resolution	0.05°C
Precision	± (1%+2 LSB) of the value
Response time at 90%	≤ 10 s

Tooxy  
Fast thermometer sensor  
Ref :  
482 204

## 4 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

### Free service \*

**+ 33 (0)2 32 29 40 50\*\***

\* Call cost not included

\*\* Only for call from foreign countries