LM 76

Lightmeter

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English

Thank you for purchasing this lightmeter.
To get the best service from this instrument :
• read these instructions carefully ;
• respect the safety precautions detailed.

1. Safety Precautions

- Do not operate the meter under environment with explosive or combustible gas, steam or filled with dust.
- To avoid reading incorrect data, replace the battery immediately when the symbol «  » appears on the LCD.

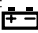
2. Features

- Overload indication : LCD will show « **OL** » in the left highest position.
- Display sampling rate : x 2.5 / sec.
- Cosine angular corrected
- Measuring light source included in the visible range
- Measuring intensities of illumination in lux or footcandles
- Length of wiring for light sensor : approx. 1.5 m
- Accessory : instruction manual, carrying case, 9V-battery

3. Description (see fig.)

1. Display (LCD) : display measurements and function symbols
2. Power button : ON/OFF switching
3. MAX button : lock up MAX or MIN, value of LCD. Press « **M-H** » again to resume normal measuring.
4. LUX/FC button : Selection of LUX and FC
5. DATA HOLD button : the reading data shown on the LCD can be locked while pressing the button.
6. RANGE button
7. Light sensor

3. Battery replacement

If the symbol «  » appears on the LCD, replace the battery immediately.

1. Remove the battery cover.
2. Replace the battery.
3. Install the battery cover.

5. Operating Instructions

1. Press the «  » button.
2. Remove the cover of the light sensor. Put the light sensor at the spot where the testing of source light is to be conducted. Auto testing will then be conducted by the meter. Read the value once stable.
3. When « **OL** » is shown at the highest position, at the left lateral side, overload is indicated. Then choose another range that is located at a comparatively higher position.
4. If you want to keep the reading value on the LCD permanently after testing, press the « **D-H** » key where by the reading value will be locked permanently up. Press the « **D-H** » key once again when you want to remove the previous locking.
5. After testing, put the cover of the light sensor back to its former position and turn off the switch.
6. After testing is completed, the indication value should be **000**, no matter what its range would be after pulling the cover of the light receiver back to its former position. Adjust **0 ADJ** to enable LCD to indicate **000**, if zero cannot be recovered. If zero still cannot be recovered, it means the meter breaks down.

Attention

- When the meter is not in use, keep the cover of the light sensor in its place to avoid the photodiode from wearing out.
- When it is not in use for a long time, take the battery away and avoid keeping it in a place of high temperature and humidity.

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