Méthode ESAO[®]

ESAO[®] Method

**Ref :
A453046**

Français – p 1

English – p 2

Version : 7011

Additif filtre antibactérien
Additional information antibacterial
filter

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Aware of the health and safety rules that are essential for biology lab work, JEULIN has included an antibacterial filter in its metabolic measurement systems.

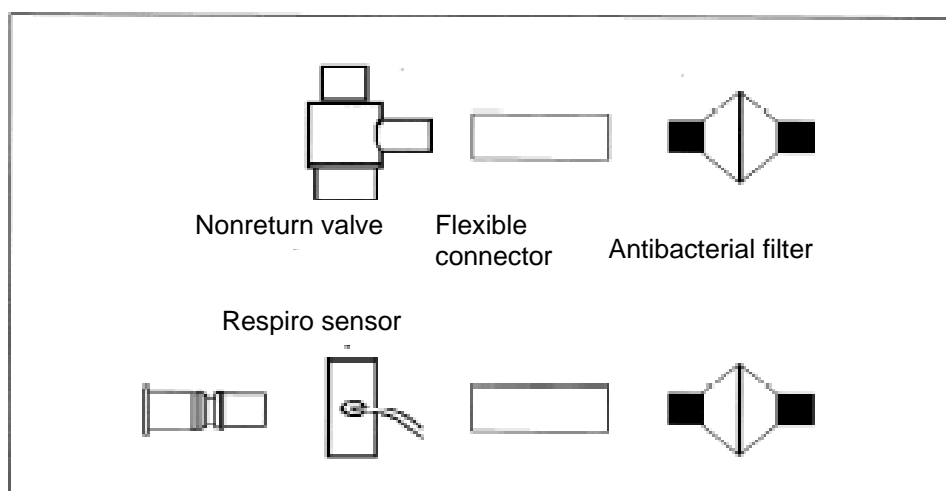
This filter achieves a major objective:

99,9 percent antibacterial and 99,7 percent antiviral protection, allowing the use in communities without the need for sterilization.

1 Installation

For RESPIHOM experiments, insert the filter between the mouthpiece and the nonreturn valve.

For RESPIRO experiments, insert the filter between the mouthpiece and the turbine

2 Precautions for use

The filters are intended for strict personal use. A method for using them may involve providing each student with a filter and identifying each filter with the student name by means of a sticker. The filters cannot be cleaned.

Decorative graphic in the top-left corner consisting of two overlapping circles, one orange and one grey.

Méthode ESAO[®]

ESAO[®] method

**Ref :
453 046**

Français – p 1

English – p 3

Version : 7011

Filtre Antibactérien
Antibacterial Filter

Decorative graphic in the bottom-right corner consisting of two overlapping circles, one orange and one grey.

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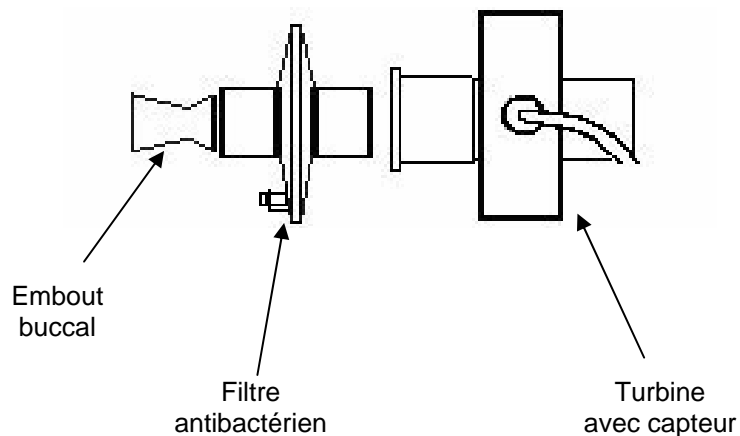
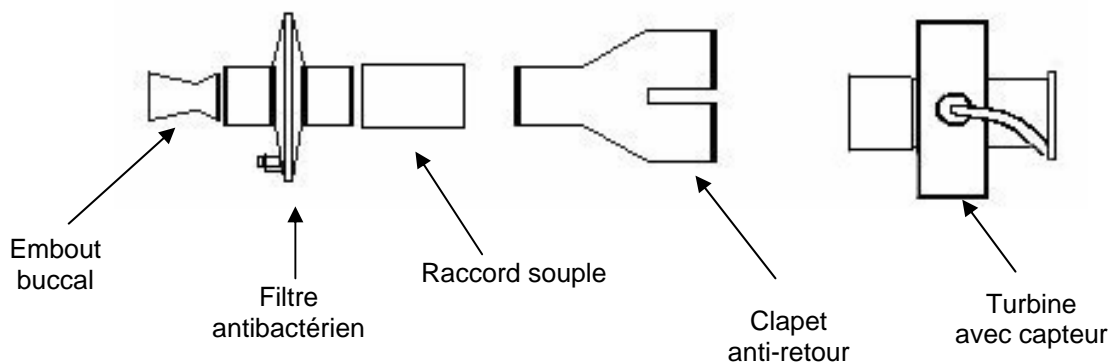
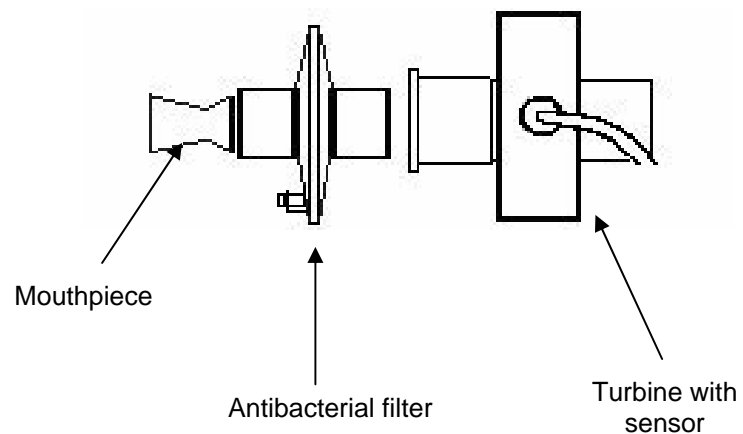
1 Description


In compliance with the health and safety rules necessary for biology TPs, this filter provides a 99.9% effective level of protection against viruses and bacteria.

2 Assembly

For the ventilation TP, insert the filter between the mouthpiece and the turbine (male side of filter = mouthpiece side).

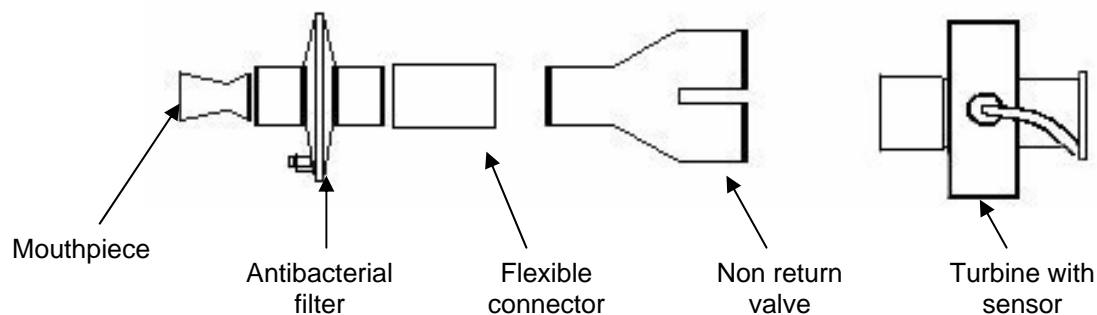For the human respiration TP, insert the filter between the mouthpiece and the non return valve (male side of filter = mouthpiece side). The non return valve is connected to the rest of the device by a flexible connector.

3 Precautions for use

Do not remove the green plug on penalty of obtaining false measurements.

The filter cannot be cleaned. Each filter is **strictly for personal use.**

4 After-Sales Service

For any additional information, please contact:

JEULIN - TECHNICAL SUPPORT
Rue Jacques Monod
BP 1900
27 019 EVREUX CEDEX FRANCE
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service * :
+ 33 (0)2 32 29 40 50

** Call cost not included*