



Test

When several Groomy® are connected to the network, one click on this button will make the selected control lamp flashing in orange.




Customize

To access this window, a password is required.

By default, the password is « expert »

But we advise you to modify it via the following button

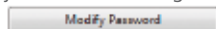


To customize your Groomy®:

- Each Groomy® device has its own name: GROOMY is the name by default.
- If several Groomy® devices are connected to the same network, customization will be necessary to be able to identify the Groomy® you want to drive and to avoid conflicts.
- Click on Name or Comment then fill in the different fields.



Settings of your Groomy®:

The Groomy® device is capable of autoconfiguration: however, the network administrators who want to set a particular IP address can do so by clicking on the following button

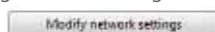

Please note:

If you forget your password or in case of settings errors of your Groomy®, the interface can be reset with the settings (name and password) by default.

To reset the interface:

Once the interface is switched off, click and hold down the button next to the power input connector 6V. Then, switch on the interface again. Hold it down about 10 sec before releasing the button.

→ Run again the application « Groomy manager »



Find all the information you need regarding your Groomy:

- Detailed characteristics
- List of compatible accessories
- List of compatible softwares
- Tips on how to use your Groomy
- Firmware updates
- Examples of some achievements



Guide for using Groomy with an Ethernet cable



- Programmable Interface
 - Standalone mode
 - Data acquisition



Throughout this guide, you will discover how simple the Groomy interface is to use.

1

Contents



- 1 Groomy interface
- 1 Power Supply 7V 2A
- 1 RJ45 connector (length 2m)
- 1 start up and maintenance manual

2

How to connect Groomy

1) Connect the plug of the processor on the Groomy device.

2) Plug in the power cord into the processor.



3) Plug in mains power plug. When this is complete, the green « status » light should flash.

4) Connect the RJ45 lead to the Groomy interface then the other end of the lead on a Ethernet network jack.

3

Installation of Groomy Manager

Groomy® interface can be connected to all kind of networks and be driven by many types of softwares
 → All devices (computer and Groomy device) must be connected to the same network.
 Groomy manager will act as a switch to assign a Groomy to a software.
 Download it on www.jeulin.com and follow the instructions.

4

Use of Groomy Manager

(Your Groomy device must be connected to the network and operated with the power supply)

In the menu Start → All the programs, → Groomy, run the Groomy manager application.
 ✂ Groomy manager

You get the following screen with the list of all Groomy devices that have been detected.



Select your Groomy interface among the displayed list.



5

The first time you connect




OK

Once your Groomy® is selected, validate to launch your application.

In our case, since we did not launch any application, the validation will execute your web browser with the display of the HTML page contained in your Groomy®
 → All the inputs/outputs can thus be driven manually.
 If you have Groomy® sensors/actuators, connect them and test them on your Groomy®.

You get then the following screen:

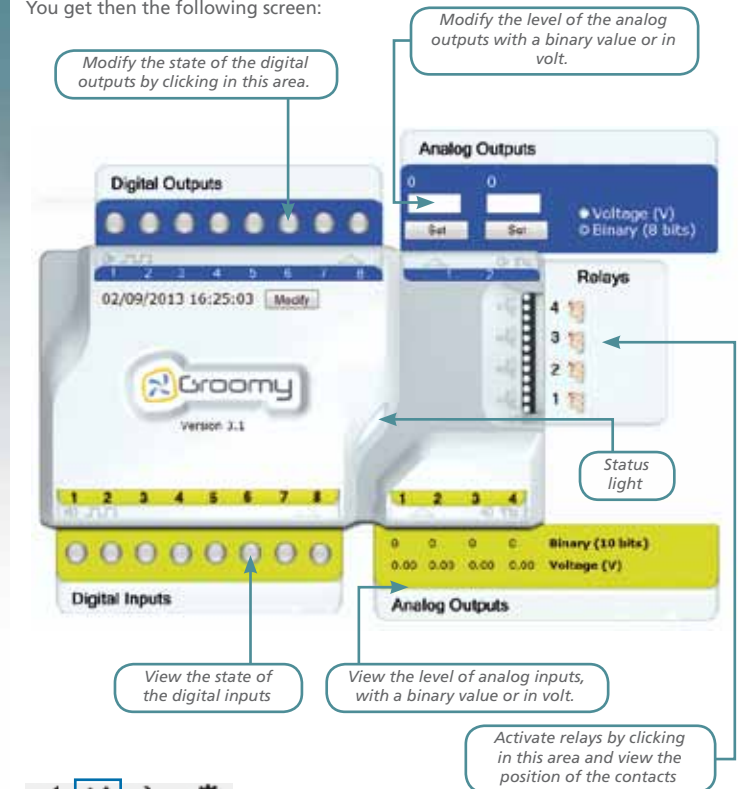

Cancel

→ Close Groomy® manager