Méthode ESAO[®]

ESAO[®] method

**Réf :
472 068**

Français – p 1

English – p 3

Version : 0111

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TRMS voltmeter sensor

1 Généralités

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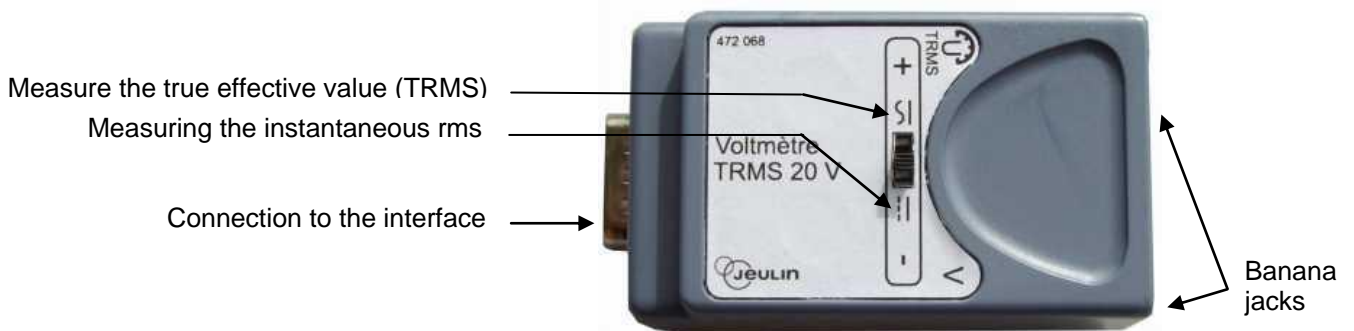
1 General

1.1 Presentation

This sensor measures the instantaneous and rms value of a voltage up to 20 V with the VTT, PRIMO and Syneo consoles. Measurements are made through a differential input with safety banana plugs.

1.2 Recognition by the VTT console

You must update your system before you can use recognition by the VTT console. You can make this update on the www.jeulin.fr internet site or using the CDROM delivered with the sensor.

This update is done connected to the PC using the Espace-VTT software:

- Connect the console to your PC
- Open the VTT space by clicking on the  icon present on your workstation
- Click on Install 
- Install the Biosflash software.

It can then be used with VTT:

- In standalone mode
- Or
- Connected to a computer (you will then need scientific workshop acquisition software).

1.3 Recognition by the PRIMO console

On the PRIMO console, the scientific workshop general-purpose software version 2.3 or later recognises the TRMS voltmeter sensor.

1.4 Recognition by the SYNEO console

The basic version of your SYNEO console recognises the TRMS Voltmeter sensor.

2 Technical characteristics

Instantaneous measurement range	From -20 V to +20 V
Effective measurement range	From 0 V to +20 V
Instantaneous measurement resolution	10 mV
Effective measurement resolution	5 mV
Instantaneous measurement precision	1 % \pm 2 digits
Effective measurement precision	2 % \pm 3 digits
Instantaneous measurement passband	From 0 to 8 kHz
Effective measurement passband	From 85 Hz to 10 kHz
Input impedance	More than 10 M Ω
Protection	Up to 40 V
Common mode voltage	40 V Max

3 After-Sales Service

For any repairs, adjustments or spare parts, please contact:

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

Free service *

+ 33 (0)2 32 29 40 50**

* Call cost not included

** Only for call from foreign countries