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## *Test and Measurement*

Réf :  
281 486

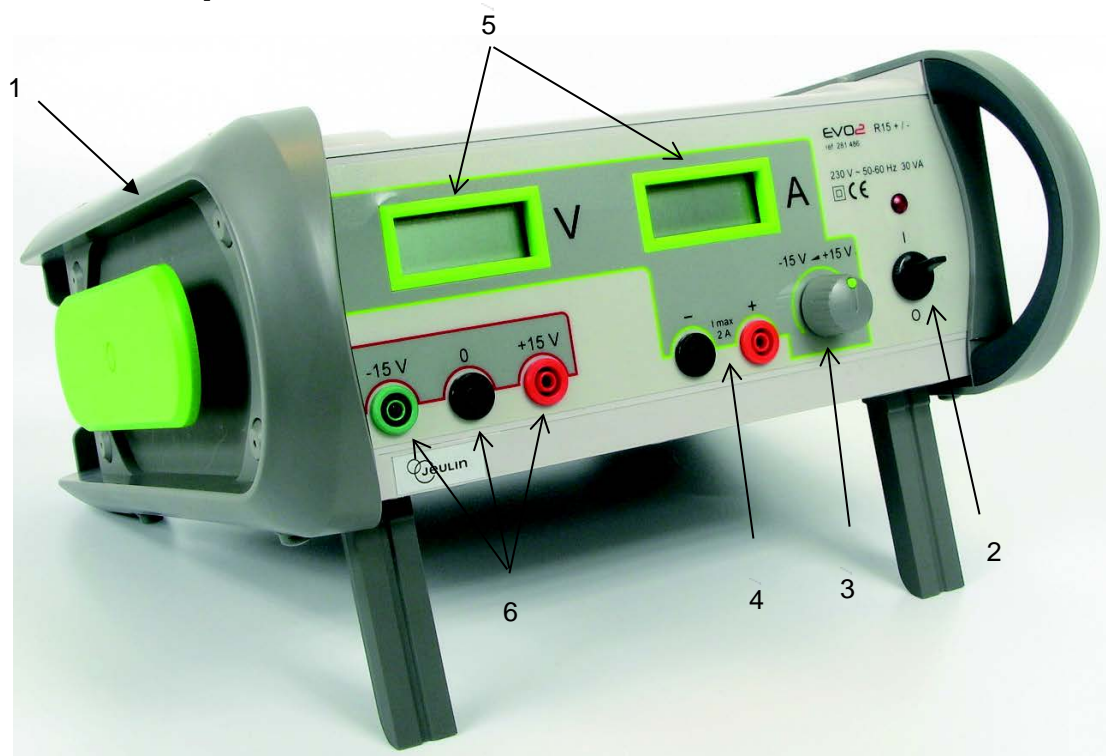
Français – p 1

English – p 3

Version : 6102

**Alimentation EVO2<sup>®</sup> R15 $\pm$  / 2 A**  
***EVO2<sup>®</sup> R15 $\pm$  / 2 A power supply***

## 1 Description

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**0 825 563 563\***

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## 1 Description

- 1 – Plastic box with double insulation
- 2 – On / off switch
- 3 – Voltage adjustment -15 V to +15 V
- 4 – Output sockets -15 V to +15 V
- 5 – External control switch
- 6 – External input sockets

## 2 Setting up

- Connect the power supply to the 230 V mains supply.
- Set the voltage and current buttons to the left (minimum value).
- Connect your circuit (lamp, motor, voltage divider, etc.) to the selected output sockets.
- Toggle the master switch (2) to " I ". The power-on indicator lamp and the 2 displays light on.
- Adjust voltage and current to desired values.

## 3 Precautions for use

- Avoid using the power supply in direct sunlight or close to a heat source.
- Never block the ventilation vents when the power supply is operating.
- If this device is used in a manner not mentioned in these instructions, the safety ensured by the device may be compromised.

## 4 Technical characteristic

- **Adjustable output** from -15 V to +15 V / 2 A,
- **Fixed output** -15 V and +15 V / 1 A,
  - Voltage drop less than 20 mV,
  - Residual ripple less than 10 mV.

**Note** that the GND of the 2 fixed outputs and the GND of the adjustable output (2 black sockets) are electrically separated. They can be linked with a banana lead.

- 2 LCD displays 2.000 points indicate output voltage and current,
- Mains voltage: 230 V  $\pm 10$  %, 50-60 Hz,
- Class II double isolation (no ground connexion),
- Consumption: 150 VA maxi,
- Primary protection with 1 A fuse slow blow (not reachable for user),
- Housing dimension: 305 x 305 x 125 mm,
- Weight: 4.5 kg.

This device complies with:

- The Low Voltage directive 2006/95/CE ; applied standard : EN 61010-1:2010,
- The EMC directive 2004/108/CEE ; applied standards : EN 60326-1.

### 4.1 Environmental conditions for use of the device

- For indoor use only,
- Altitude less than 2000 meters,
- Operating temperature from 15 °C to 40 °C,
- Maximum relative humidity of 80 % up to 31 °C and 50 % up to 40 °C,
- Mains supply variations of  $\pm 10$  %,
- Installation category II, pollution level 2.

### 4.2 Care and maintenance

In case of a malfunction:

- Check the presence of a mains voltage (230 V), the red indicator lamp under the on / off switch is on.
- Check your set-up so as to detect any possible short circuits.

If the device still does not work, contact the After Sales Service.

In any cases, **do not try to open the device.**

Store the original packaging; it can be used to bring the device to our workshops for maintenance, if necessary.

## 5 After-Sales Service

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

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# Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediately to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

## Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.  
\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux