**Ex.A.O.**

Capteur Primo / VTT

***Datalogging***

*Primo / VTT sensor*

Réf :  
472 010

Français – p 1

English – p 3

Version : 8109

**Capteur pression 0/2000 hPa**

***Pressure sensor 0/2000 hPa***

## 1. Description

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**0 825 563 563\***

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## 1. Description

This sensor enables measurement of absolute pressure from 0 to 2000 hPa. It is supplied with a 1 m length of tube.

Measurements are made by means of a semiconductor-based sensor.  
This sensor is used with the VTT® and Primo® interfaces.

## 2. Use

### 2.1 Connecting the sensor to the console

- Introduce the sensor aligned with the measurement channel chosen.
- Locate the sensor in the console's guide slots.
- Follow the console's guide slot system.
- Locate fully home.

### 2.2 Use

- Connect the sensor.
- Switch on the console.

## 3. Technical features

<b>Measurement range</b>	From 0 to 2000 hPa
<b>Resolution</b>	2 hPa
<b>Accuracy</b>	2% $\pm$ 4 hPa
<b>Environnement</b>	Non-corrosive gases (not to be used in liquids)
<b>Maximum pressure</b>	4000 hPa
<b>Recognition code</b>	8

## 4. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

**JEULIN – TECHNICAL SUPPORT**  
468 rue Jacques Monod  
CS 21900  
27019 EVREUX CEDEX FRANCE

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# Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediatly to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

## Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.

\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux