**Ex.A.O.**

Capteur Foxy / Tooxy / AirNeXT

***Data logging***

*Sensor*

Réf :  
482 076

Français – p 1

English – p 5

Version : 8106

**Capteur CO<sub>2</sub> Air**  
***Air CO<sub>2</sub> Sensor***

## 1. Description

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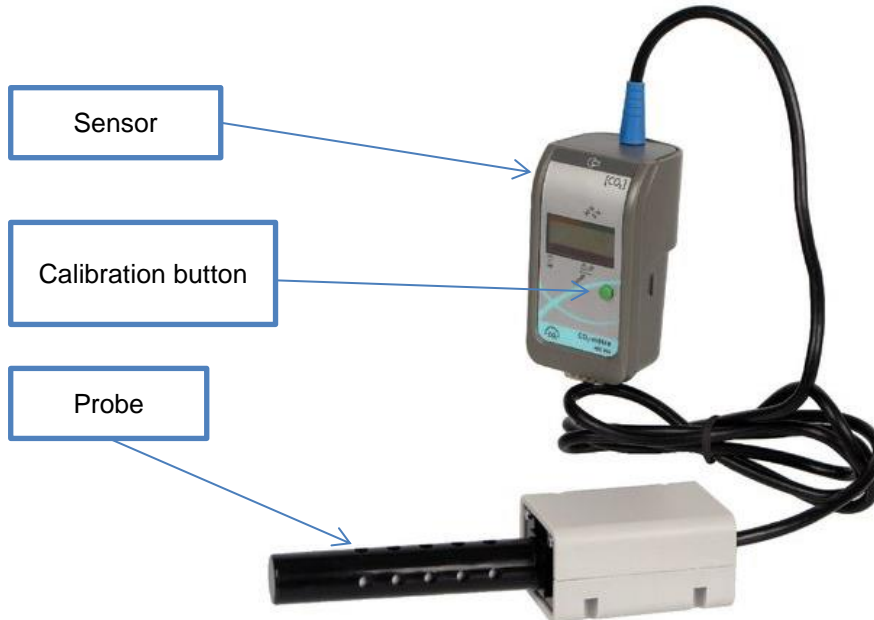
**0 825 563 563\***

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## 1. Description

The Air CO<sub>2</sub> sensor measures CO<sub>2</sub> concentration in the air. It can be used for data capture or a direct reading via its display.

## 2. Use

### 2.1 Warning

Do not immerse the probe in liquid or it may be damaged permanently.

**For a better precision, switch on the probe 5 minutes before any manipulation.**

### 2.2 Interface

This sensor is used with Foxy®, AirNeXT® and Tooxy® consoles.

### 2.3 Inserting the sensor

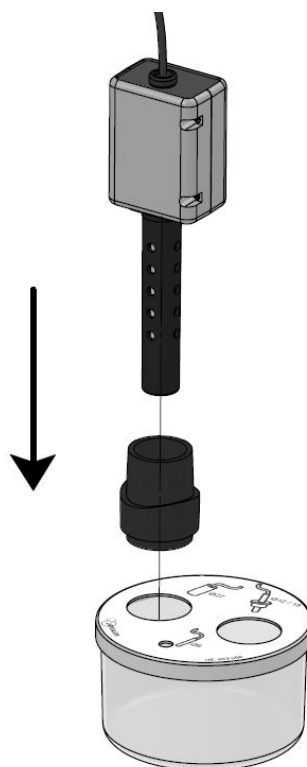
- Place the sensor facing the console connector
- Slide the sensor following the console guide
- Push until it clicks

### 2.4 Inserting the probe


The sensor is provided with a probe adapter. It allows the probe to be used with the following products

- Animal respiration device (ref: 453 066)
- Human respiration device (ref: 453 143)
- Human metabolism device (ref: 453 144)

For the animal respiration device:

For human respiration devices
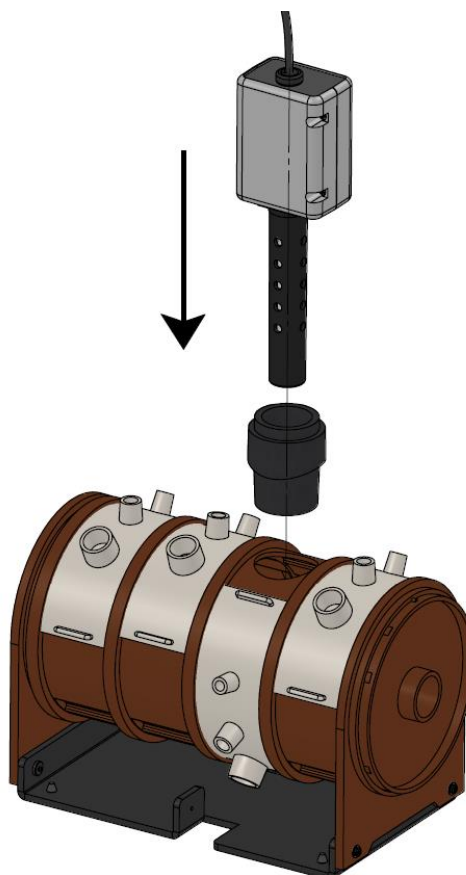
## 2.5 Use in measurement mode

- The sensor has a 1 minute warm-up period.
- During this time, the display is inactive. Once the warm-up period is complete the CO<sub>2</sub> rate appears on the screen.

### 0.04% adjustment:

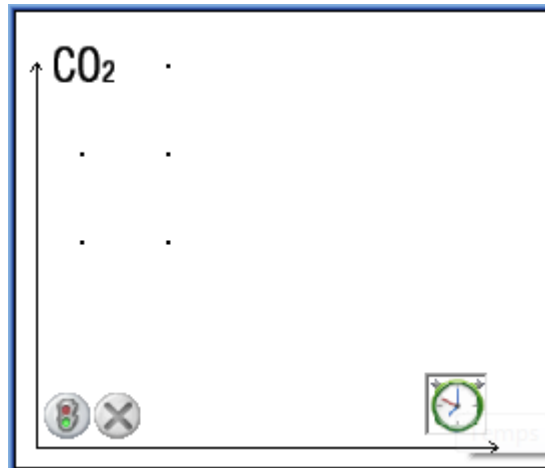
- Push and hold the calibration button
- Wait until the value appears on the screen
- Push the button briefly to confirm the calibration

**Note:** it's possible that the posted value differs slightly from 0.04 %. In this case, begin again the procedure of calibration.

## 2.6 Use with a console

The buttons are inactive when the sensor is connected to a console. Adjustments are made via the software.

The CO<sub>2</sub> sensor is represented by the **CO<sub>2</sub>** icon in the Data Science Studio general module. To complete a capture, drag the icon along one of the axes. For example, to capture the CO<sub>2</sub> rate based on time, drag the CO<sub>2</sub> icon along the Y axis and the time icon along the X axis.

### 0.04% adjustment:

- Place the sensor icon on the Y axis
- Click on the "Adjustment" tab
- Click on Start
- Enter 0.04% in the box
- Wait until the red rectangle turns green.
- Click on "Adjust"

### 3. Technical characteristics

Warm-up period	30s - 1 minute
Measurement range	0 - 10% of CO <sub>2</sub>
Measurement accuracy	at 10% of CO <sub>2</sub> : $\pm 0,8\%$ of CO <sub>2</sub> at 5% of CO <sub>2</sub> : $\pm 0,03\%$ of CO <sub>2</sub> at 0,2% of CO <sub>2</sub> : $\pm 0,01\%$ of CO <sub>2</sub> at 0,1% of CO <sub>2</sub> : $\pm 0,006\%$ of CO <sub>2</sub>
Probe useful life	>5 years

### 4. After-sales service

The device is under a 2-year guarantee, it must be sent back to our workshops.  
For any repairs, adjustments or spare parts please contact:

**JEULIN – TECHNICAL SUPPORT**  
468 rue Jacques Monod  
CS 21900  
27019 EVREUX CEDEX FRANCE

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## Direct connection for technical support

A team of experts  
at your disposal  
from Monday to Friday  
(opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request  
immediately to provide you  
with the right answers regarding  
your activity field : Biology, Physics,  
Chemistry, Technology.

### Free service\*

**+33 2 32 29 40 50\*\***

\* Call cost not included.  
\*\* Only for call from foreign countries.

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SAS au capital de 1 000 000 € - TVA intracommunautaire FR47 344 652 490 - Siren 344 652 490 RCS Evreux